Ultramikrotome (PowerTomes, RMC)

Ultramicrotomes (PowerTomes, RMC)

Kryokammer (LN Ultra, RMC)

Cryo Chamber (LN Ultra, RMC)

Array Tomographie (ATUMtome, RMC)

Array Tomography (ATUMTOME, RMC)

RMC Boeckeler PowerTome Serie

RMC Boeckeler PowerTome series

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The RMC PowerTome series is a highly stable and robust ultramicrotome for biological, material science and clinical applications. The PowerTome is also the basis for the ATUMtome, a section collection system for large volume 3D reconstructions. Therefore, the components are built and tested for high work load and long-term stability. The PowerTome is available in various upgradeable models.

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The PT XL is the basic model with tactile control unit for the ultramicrotome operation and is eminently suitable for routine work. This model is upgradeable to all versions including the LN Ultra attachment for cryosectioning and the ATUMtome for automated serial section collection.

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All Powertomes are capable of accommodating glass knives up to 12mm wide, triangular tungsten carbide knives and diamond knives of any commercial brand.

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► Power Drive

Motorized cutting stroke, supports a homogeneous sectioning without vibrations even for large, hard and inhomogeneous specimen blocks.

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► Intergrated Handrest

For resting the hands, e.g. during manipulation of ultra-thin sections supporting an ergonomic work environment (removable).

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► **Cutting Speed**

0.1 to 100mm/s cutting speed adjustable in 0.1mm/s steps throughout whole speed range.

► **Scan and Tilt**

A small knob on the microscope turret enabling to scan the knife edge area without turning the entire microscope even at high magnifications.

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PT PC with HD video package and view for training and documentation

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The PT PC has all the features of the PT XL and is further equipped with PC control and camera for live image acquisition and a video and snapshot function. The PC control enables documentation and archiving of user settings and protocols. The online support using internet capabilities is included, for a direct link to specialists to assist troubleshooting.

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The high definition video package to monitor the sectioning process in real-time is displayed at the center of the PC monitor. An image capture system allows videos or still photographs to be taken at any time for data sharing in

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the lab or for collaboration as well as for publications or archiving.

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The ultramicrotome can be controlled with the touch screen monitor the digital tactile controller or with both.

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► **Dual Drive**

Touch screen PC and digital tactile controller.

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► **Video Monitoring**

The center of the touch screen displays the sectioning process in real time. Essential for user trainings and demonstrations. With recording function for acquisition of video and snap shots for documentation.

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► **VisuTrac**

Define and adjust cutting window during sectioning with a single mouse click.

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► **Digital Measuring Tool**

Define and store sample and block dimensions.

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► **Trimming Function**

Automated trimming into the sample block with defined depth (max. 200µm).

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► **Integrated Report Generator**

Direct read-out on PC to user storage via ethernet, without the need of portable data storage (e.g. USB stick).

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► **Online Trouble Shooting**

Skype access to RMC Service engineers.

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
PT 3D - first 1mm advance ultramicrotome

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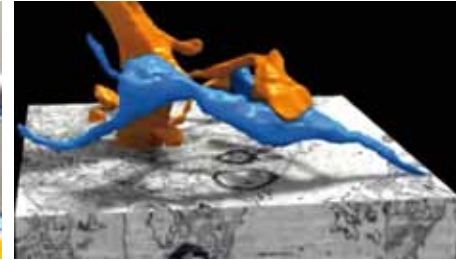For decades, a specimen advance of max. 200µm was the standard in ultramicrotomy. With the newly designed cutting mechanism, the PT 3D sets a new standard and is the first ultramicrotome with the ability to cut ultrathin sections up to one millimeter into a sample without resetting the advance mechanism.

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The unique 1mm specimen advance is not only beneficial for ultramicrotomy applications such as 3D reconstruction of materials, tissues, cells or cellular structures using array tomography, but also for those involving frequent trimming with the highest precision, like for the sample preparation for cryosectioning.

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Identical to the PT PC the PT 3D has a high definition video package to display and control the section process in real-time on the computer monitor. The PC software further includes all other functions, e.g. the image capture system, the digital measurement tools and more.

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► Reduced Handling

The rate of interfering with knife and sample because of resetting the advance is reduced by five times.

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► Every section is sacred

For 3D reconstruction, it is crucial to reduce the number of lost sections to a minimum. The main risk of losing a section is the re-approach to the knife edge.

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► Less Knife Damage

Most damage to a diamond knife occurs during approach of knife and sample.

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
► **RMC Specimen Holders**

Made of sturdy stainless steel with side entry for backlighting the sample in the block. Backlighting also possible for pre-trimming with the trimming block. With diameter of 10 mm, compatible with ultramicrotomes of common manufacturers.

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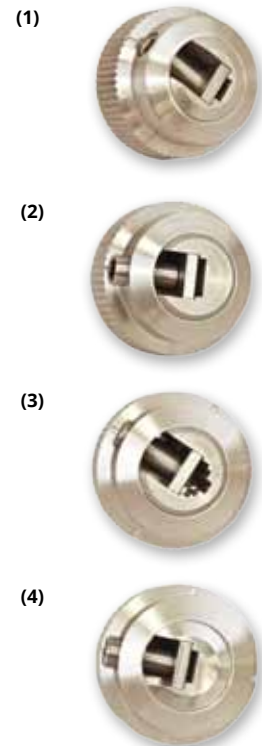► **DiATOME ultra 45°, 3mm**

Standard knife for routine work. High knife stability with higher compression due to 45° angle.

DU4530	DiATOME ultra 45° 3,0mm <i>DiATOME ultra 45° 3.0mm</i>
--------	---

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► **DiATOME ultra 35°, 3mm**

Delicate knife for thinner sections and series. Smaller compression with more fragile cutting edge.

DU4530	DiATOME ultra 45° 3,0mm <i>DiATOME ultra 45° 3.0mm</i>
--------	---

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► **Vibration Isolation table by TMC**

Compact design and optimized ergonomics, compressed air connection required.

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LN Ultra - Cryo chamber for highest temperature stability

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Sectioning of both biological and industrial materials at cryogenic temperatures can be achieved with the LNUltra a liquid nitrogen-cooled cryo attachment for the PowerTome.

Capable of holding temperatures down to -180°C , the stable environment of the LNUltra is perfect for sectioning and planing a variety of samples including rubber and soft polymers, as well as biological tissues and cells, e.g., for immunolabelling using the Tokuyasu technique or CEMOVIS (Cryo Electron Microscopy of Vitreous Sections).

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Due to the low Liquid Nitrogen (LN2) consumption, a small 12-liter table-top Dewar is sufficient to provide a stable sectioning environment for around 6 hours, while saving LN2 costs.

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In combination with the PowerTome 3D, sample trimming and cutting can be achieved without resetting the specimen advance and without the need for realigning sample and knife.

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► Low LN2 Consumption

Approx. 2L/h with use of a space-saving 12L tabletop dewar.

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► High Temperature Stability

From $+20^{\circ}\text{C}$ to -180°C due to use of exclusive materials.

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► In-Chamber Specimen Freezing

Open space in the chamber for easy freezing or mounting the specimen on the pin.

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► **Easy Chamber Access**

Reduced depth of chamber facilitates sample handling and manipulation.

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► **Workstation**

Lay up tools and hands for working at the chamber without becoming cold.

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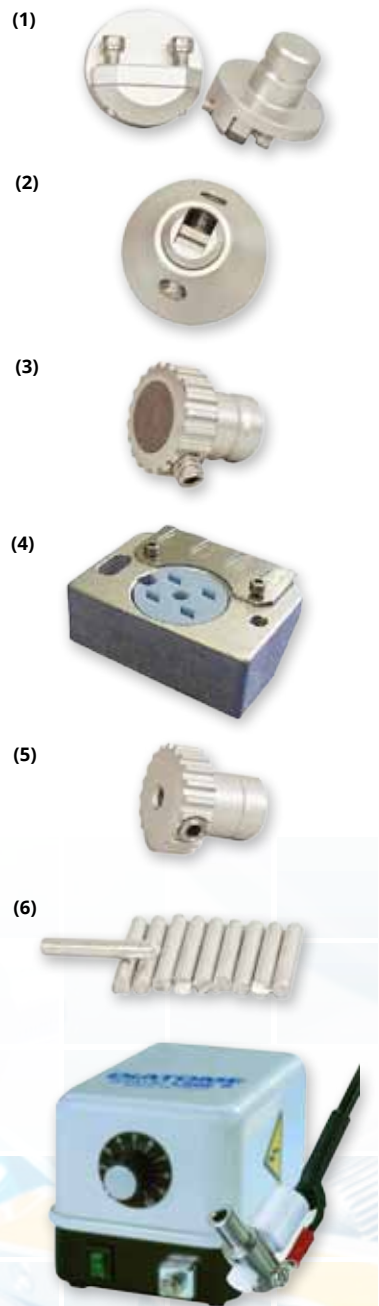
Reduced electrostatic charging of specimens

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DCIMM3530

► **DiATOME cryo immuno 35°, 3mm for Tokuyasu**

Specially manufactured for sample preparation using the Tokuyasu technique

DiATOME cryo immuno 35° 3,0mm, für Tokuyasu

DiATOME cryo immuno 35° 3,0mm, for Tokuyasu

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DCO3530

► **DiATOME cryo dry 35°, 3mm**

for dry sectioning of material samples, e.g. polymers and films

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Diatome trim 45°, for precise trimming of sample blocks at room and cryo temperatures

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E71952-01

► **Blades**

*Rough pre-trimming of sample blocks. **Caution: Only use cleaned blades***

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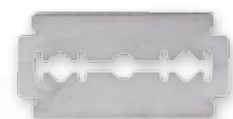
Single Edged Razor Blades for Clean Rooms, Surgical Stainless Steel

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Double Edged Razor Blades, Feather®, Carbon Steel

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Dumont Style 5, straight Tweezer for handling of TEM-Grids and samples

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Dumont Style 7 curved Tweezer for handling of TEM-Grids and samples

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Tech@ScienceServices.de

Any further Questions?

If you have any further questions regarding a product or the field of microtomy, please do not hesitate to contact us. We also offer on-site demonstrations.

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Science Services GmbH